Code No. RR916

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Takara

Real Time PCR
Chicken DNA Detection

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Thermal Cycler Dice™ Real Time System III (Code No. TP950/TP970/TP980/TP990)

Thermal Cycler Dice Real Time System *Lite* (Code No. TP700/TP760)

Applied Biosystems 7300/7500/7500 Fast Real-Time PCR System and StepOnePlus Real-Time PCR System (Thermo Fisher Scientific)

LightCycler 96/480 System (Roche Diagnostics)

CFX96 Real-Time PCR Detection System (Bio-Rad)

Smart Cycler System/Smart Cycler II System (Cepheid)

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	Target	Reporter	Quencher
Probe Mix	Chicken	FAM	Eclipse
	Internal Control	HEX	Eclipse

● Control DNA for Chicken 及 Internal Control DNA 扩增片段大小

DNA 种类	扩增片段大小
Control DNA for Chicken	114 bp
Internal Control DNA	92 bp

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1. Real Time PCR Bovine DNA Detection Kit (Code No. RR910)
2. Real Time PCR Ovine DNA Detection Kit (Code No. RR911)
3. Real Time PCR Porcine DNA Detection Kit (Code No. RR912)
4. Real Time PCR Bovine and Ovine DNA Detection Kit (Code No. RR913)
5. Real Time PCR Rabbit DNA Detection Kit (Code No. RR914)
6. Real Time PCR Mammalian DNA Detection Kit (Code No. RR915)
7. Real Time PCR Cat DNA Detection Kit (Code No. RR926)
8. Real Time PCR Martes DNA Detection Kit (Code No. RR928)
9. Real Time PCR 18s DNA Detection Kit (Code No. RR930)
10. Real Time PCR Bos grunniens DNA Detection Kit (Code No. RR931)
11. Real Time PCR Bos taurus DNA Detection Kit (Code No. RR932)
12. Real Time PCR Bubalus bubalis DNA Detection Kit (Code No. RR933)
13. Real Time PCR Duck DNA Detection Kit (Code No. RR934)
14. TaKaRa MiniBEST Universal Genomic DNA Extraction Kit Ver.5.0 (Code No. 9765)
15. DNA Isolation Reagent for Meat and Meat Products (Code No. 9178)

TaKaRa Ex Taq is a registered trademark of Takara Bio Inc.

Thermal Cycler Dice is a trademark of Takara Bio Inc.

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